

miRCURY LNA miRNA miRNome PCR Panels

Human Panel 6, V6 Product Data Sheet

Cat. no. 339322 YAHS-306X

For mature miRNA expression profiling using real-time PCR

Format	Suitable real-time cyclers	Plate	Cat. no.
E	Applied Biosystems® models 7900HT (384-well block), ViiA™ 7 (384-well block); Bio-Rad CFX384™	384-well	YAHS-306XE
G	Roche® LightCycler® 480 (384-well block)	384-well	YAHS-306XG

Description

The miRCURY LNA miRNA miRNome PCR Panels profile the expression of hundreds of the most abundantly expressed and best characterized miRNA sequences in the a given species-specific miRNome as annotated in miRBase (www.mirbase.org). Although they are well characterized, each of these miRNA sequences can regulate one to several messenger RNA transcripts, and conversely one mRNA can be regulated by one to several miRNA sequences. Therefore, the complex role that any given known miRNA sequence plays has yet to be completely defined. Use of the miRCURY LNA miRNA miRNome PCR Panels maximizes the likelihood of discovering miRNA sequences whose expression patterns correlate with the biological phenotypes under study. A set of controls included on each plate enables data analysis using the $\Delta\Delta C_q$ method of relative quantification, assessment of reverse transcription performance, and assessment of PCR performance. The miRCURY LNA miRNA miRNome PCR Panels enable easy and reliable, SYBR Green-based real-time PCR analysis of the expression of miRNA sequences and discovery of those most important to the research area of interest.

For further details, consult the *miRCURY LNA™ miRCURY LNA™ miRNA Focus PCR Panels Handbook*.

Array Layout

15	16	17	18	19	20	21	22	23	24
hsa-miR-1250-3p	hsa-miR-1251-3p	hsa-miR-1252-3p	hsa-miR-1266-3p	hsa-miR-1273n-3p	hsa-miR-128-1-5p	hsa-miR-128-2-5p	hsa-miR-1288-5p	hsa-miR-1296-3p	hsa-miR-1298-3p
hsa-miR-216b-3p	hsa-miR-2276-5p	hsa-miR-3151-3p	hsa-miR-3192-3p	hsa-miR-328-5p	hsa-miR-329-5p	hsa-miR-370-5p	hsa-miR-372-5p	hsa-miR-383-3p	hsa-miR-3912-5p
hsa-miR-5189-3p	hsa-miR-520g-5p	hsa-miR-548ba	hsa-miR-548e-5p	hsa-miR-548f-5p	hsa-miR-548i-3p	hsa-miR-5699-5p	hsa-miR-579-5p	hsa-miR-580-5p	hsa-miR-585-5p
hsa-miR-6728-3p	hsa-miR-6730-5p	hsa-miR-6731-3p	hsa-miR-6732-3p	hsa-miR-6732-5p	hsa-miR-6733-3p	hsa-miR-6733-5p	hsa-miR-6734-3p	hsa-miR-6734-5p	hsa-miR-6735-5p
hsa-miR-6744-3p	hsa-miR-6744-5p	hsa-miR-6746-3p	hsa-miR-6747-3p	hsa-miR-6748-3p	hsa-miR-6748-5p	hsa-miR-6749-5p	hsa-miR-6750-5p	hsa-miR-6751-3p	hsa-miR-6751-5p
hsa-miR-6762-5p	hsa-miR-6763-5p	hsa-miR-6764-3p	hsa-miR-6764-5p	hsa-miR-6765-5p	hsa-miR-6766-5p	hsa-miR-6767-3p	hsa-miR-6767-5p	hsa-miR-6768-3p	hsa-miR-6768-5p
hsa-miR-6775-5p	hsa-miR-6776-3p	hsa-miR-6776-5p	hsa-miR-6777-5p	hsa-miR-6778-3p	hsa-miR-6778-5p	hsa-miR-6779-3p	hsa-miR-6779-5p	hsa-miR-6780a-3p	hsa-miR-6780a-5p
hsa-miR-6790-3p	hsa-miR-6790-5p	hsa-miR-6791-3p	hsa-miR-6792-3p	hsa-miR-6792-5p	hsa-miR-6793-5p	hsa-miR-6794-3p	hsa-miR-6794-5p	hsa-miR-6795-3p	hsa-miR-6795-5p
hsa-miR-6805-5p	hsa-miR-6806-3p	hsa-miR-6806-5p	hsa-miR-6807-3p	hsa-miR-6807-5p	hsa-miR-6808-3p	hsa-miR-6808-5p	hsa-miR-6809-3p	hsa-miR-6809-5p	hsa-miR-6810-5p
hsa-miR-6820-5p	hsa-miR-6821-3p	hsa-miR-6821-5p	hsa-miR-6822-3p	hsa-miR-6822-5p	hsa-miR-6823-3p	hsa-miR-6823-5p	hsa-miR-6824-3p	hsa-miR-6824-5p	hsa-miR-6825-3p
hsa-miR-6832-5p	hsa-miR-6833-3p	hsa-miR-6833-5p	hsa-miR-6834-3p	hsa-miR-6834-5p	hsa-miR-6835-3p	hsa-miR-6835-5p	hsa-miR-6836-3p	hsa-miR-6836-5p	hsa-miR-6837-3p
hsa-miR-6846-3p	hsa-miR-6846-5p	hsa-miR-6847-3p	hsa-miR-6847-5p	hsa-miR-6848-3p	hsa-miR-6848-5p	hsa-miR-6849-3p	hsa-miR-6849-5p	hsa-miR-6850-3p	hsa-miR-6851-3p
hsa-miR-6860	hsa-miR-6861-5p	hsa-miR-6862-3p	hsa-miR-6862-5p	hsa-miR-6863	hsa-miR-6864-3p	hsa-miR-6864-5p	hsa-miR-6865-3p	hsa-miR-6865-5p	hsa-miR-6866-3p
hsa-miR-6875-3p	hsa-miR-6875-5p	hsa-miR-6876-3p	hsa-miR-6876-5p	hsa-miR-6877-5p	hsa-miR-6878-3p	hsa-miR-6878-5p	hsa-miR-6879-5p	hsa-miR-6880-5p	hsa-miR-6881-3p
hsa-miR-6890-5p	hsa-miR-6891-3p	hsa-miR-6891-5p	hsa-miR-6892-3p	hsa-miR-6892-5p	hsa-miR-6893-3p	hsa-miR-6893-5p	hsa-miR-7107-3p	hsa-miR-7108-3p	hsa-miR-7109-5p
SNORD49A (hsa)	U6 snRNA (v2)	cel-miR-39-3p	Unisp2	Unisp4	Unisp5	Unisp6	Unisp3	Unisp3	Unisp3

5	6	7	8	9	10	11	12	13	14
hsa-miR-6722-3p	hsa-miR-6722-5p	hsa-miR-758-5p	hsa-miR-892c-3p	hsa-miR-892c-5p	hsa-miR-937-5p	hsa-miR-98-3p	hsa-let-7c-3p	hsa-miR-1180-5p	hsa-miR-1199-3p
hsa-miR-1468-3p	hsa-miR-152-5p	hsa-miR-1537-5p	hsa-miR-181d-3p	gga-miR-190a-3p	hsa-miR-1910-3p	hsa-miR-208a-5p	hsa-miR-208b-5p	hsa-miR-210-5p	hsa-miR-215-3p
hsa-miR-4433b-3p	hsa-miR-4433b-5p	hsa-miR-487a-5p	hsa-miR-487b-5p	hsa-miR-489-5p	hsa-miR-494-5p	hsa-miR-500b-3p	hsa-miR-6889-3p	hsa-miR-511-3p	hsa-miR-513b-3p
hsa-miR-6516-3p	hsa-miR-6516-5p	hsa-miR-653-3p	hsa-miR-655-5p	hsa-miR-656-5p	hsa-miR-668-5p	hsa-miR-670-3p	hsa-miR-6889-5p	hsa-miR-6726-3p	hsa-miR-6726-5p
hsa-miR-6738-5p	hsa-miR-6739-3p	hsa-miR-6739-5p	hsa-miR-6740-3p	hsa-miR-6740-5p	hsa-miR-6741-3p	hsa-miR-6741-5p	hsa-miR-6742-3p	hsa-miR-6742-5p	hsa-miR-6743-5p
hsa-miR-6755-3p	hsa-miR-6755-5p	hsa-miR-6756-5p	hsa-miR-6757-3p	hsa-miR-6757-5p	hsa-miR-6758-3p	hsa-miR-6759-3p	hsa-miR-6759-5p	hsa-miR-6761-3p	hsa-miR-6762-3p
hsa-miR-6770-5p	hsa-miR-6771-3p	hsa-miR-6771-5p	hsa-miR-6772-3p	hsa-miR-6772-5p	hsa-miR-6773-3p	hsa-miR-6773-5p	hsa-miR-6774-3p	hsa-miR-6774-5p	hsa-miR-6775-3p
hsa-miR-6783-5p	hsa-miR-6785-3p	hsa-miR-6785-5p	hsa-miR-6786-3p	hsa-miR-6786-5p	hsa-miR-6787-3p	hsa-miR-6787-5p	hsa-miR-6788-3p	hsa-miR-6788-5p	hsa-miR-6789-5p
hsa-miR-6799-5p	hsa-miR-6800-3p	hsa-miR-6800-5p	hsa-miR-6801-3p	hsa-miR-6801-5p	hsa-miR-6802-3p	hsa-miR-6802-5p	hsa-miR-6803-3p	hsa-miR-6804-3p	hsa-miR-6804-5p
hsa-miR-6814-5p	hsa-miR-6815-3p	hsa-miR-6815-5p	hsa-miR-6816-3p	hsa-miR-6816-5p	hsa-miR-6817-5p	hsa-miR-6818-3p	hsa-miR-6818-5p	hsa-miR-6819-5p	hsa-miR-6820-3p
hsa-miR-6827-5p	hsa-miR-6828-3p	hsa-miR-6828-5p	hsa-miR-6829-3p	hsa-miR-6829-5p	hsa-miR-6830-3p	hsa-miR-6830-5p	hsa-miR-6831-3p	hsa-miR-6831-5p	hsa-miR-6832-3p
hsa-miR-6839-5p	hsa-miR-6840-3p	hsa-miR-6840-5p	hsa-miR-6841-3p	hsa-miR-6841-5p	hsa-miR-6842-3p	hsa-miR-6842-5p	hsa-miR-6843-3p	hsa-miR-6890-3p	hsa-miR-6845-5p
hsa-miR-6853-5p	hsa-miR-6854-3p	hsa-miR-6854-5p	hsa-miR-6855-3p	hsa-miR-6855-5p	hsa-miR-6856-3p	hsa-miR-6857-3p	hsa-miR-6857-5p	hsa-miR-6858-5p	hsa-miR-6859-5p
hsa-miR-6868-5p	hsa-miR-6869-3p	hsa-miR-6870-3p	hsa-miR-6871-3p	hsa-miR-6871-5p	hsa-miR-6872-5p	hsa-miR-6873-3p	hsa-miR-6873-5p	hsa-miR-6874-3p	hsa-miR-6874-5p
hsa-miR-6883-5p	hsa-miR-6884-3p	hsa-miR-6884-5p	hsa-miR-6885-3p	hsa-miR-6885-5p	hsa-miR-6886-5p	hsa-miR-6887-3p	hsa-miR-6887-5p	hsa-miR-6888-3p	hsa-miR-6888-5p
hsa-miR-7113-3p	hsa-miR-7113-5p	hsa-miR-7114-3p	hsa-miR-7114-5p	hsa-miR-7150	hsa-miR-7151-3p	hsa-miR-7151-5p	BlankSpot	SNORD44 (hsa)	SNORD38B (hsa)

	1	2	3	4
A	hsa-miR-6718-5p	hsa-miR-6719-3p	hsa-miR-6720-3p	hsa-miR-6721-5p
B	hsa-miR-1301-5p	hsa-miR-133a-5p	hsa-miR-1343-5p	hsa-miR-1343-p
C	hsa-miR-3928-5p	hsa-miR-410-5p	hsa-miR-412-5p	hsa-miR-433-5p
D	hsa-miR-598-5p	hsa-miR-605-3p	hsa-miR-619-5p	hsa-miR-651-3p
E	hsa-miR-6736-3p	hsa-miR-6736-5p	hsa-miR-6737-5p	hsa-miR-6738-3p
F	hsa-miR-6752-5p	hsa-miR-6753-3p	hsa-miR-6753-5p	hsa-miR-6754-5p
G	hsa-miR-6769a-5p	hsa-miR-6769b-3p	hsa-miR-6769b-5p	hsa-miR-6770-3p
H	hsa-miR-6780b-3p	hsa-miR-6781-5p	hsa-miR-6782-3p	hsa-miR-6783-3p
I	hsa-miR-6796-5p	hsa-miR-6797-5p	hsa-miR-6798-5p	hsa-miR-6799-3p
J	hsa-miR-6811-3p	hsa-miR-6811-5p	hsa-miR-6812-5p	hsa-miR-6813-5p
K	hsa-miR-6825-5p	hsa-miR-6826-3p	hsa-miR-6826-5p	hsa-miR-6827-3p
L	hsa-miR-6837-5p	hsa-miR-6838-3p	hsa-miR-6838-5p	hsa-miR-6839-3p
M	hsa-miR-6851-5p	hsa-miR-6852-3p	hsa-miR-6852-5p	hsa-miR-6853-3p
N	hsa-miR-6866-5p	hsa-miR-6867-3p	hsa-miR-6867-5p	hsa-miR-6868-3p
O	hsa-miR-6881-5p	hsa-miR-6882-3p	hsa-miR-6882-5p	hsa-miR-6883-3p
P	hsa-miR-7110-3p	hsa-miR-7111-3p	hsa-miR-7112-3p	hsa-miR-7112-5p

miRNA Table

Well	miRNA ID	Accession #	Assay Catalog #	Well	miRNA ID	Accession #	Assay Catalog #
A01	hsa-miR-6718-5p	MIMAT0025849	YP02113259	I01	hsa-miR-6796-5p	MIMAT0027492	YP02112426
A02	hsa-miR-6719-3p	MIMAT0025850	YP02111364	I02	hsa-miR-6797-5p	MIMAT0027494	YP02118253
A03	hsa-miR-6720-3p	MIMAT0025851	YP02114592	I03	hsa-miR-6798-5p	MIMAT0027496	YP02105728
A04	hsa-miR-6721-5p	MIMAT0025852	YP02110957	I04	hsa-miR-6799-3p	MIMAT0027499	YP02101545
A05	hsa-miR-6722-3p	MIMAT0025854	YP02105759	I05	hsa-miR-6799-5p	MIMAT0027498	YP02100691
A06	hsa-miR-6722-5p	MIMAT0025853	YP02113668	I06	hsa-miR-6800-3p	MIMAT0027501	YP02107834
A07	hsa-miR-758-5p	MIMAT0022929	YP02116137	I07	hsa-miR-6800-5p	MIMAT0027500	YP02112192
A08	hsa-miR-892c-3p	MIMAT0025858	YP02117971	I08	hsa-miR-6801-3p	MIMAT0027503	YP02101519
A09	hsa-miR-892c-5p	MIMAT0025857	YP02114947	I09	hsa-miR-6801-5p	MIMAT0027502	YP02115784
A10	hsa-miR-937-5p	MIMAT0022938	YP02117509	I10	hsa-miR-6802-3p	MIMAT0027505	YP02112736
A11	hsa-miR-98-3p	MIMAT0022842	YP02106353	I11	hsa-miR-6802-5p	MIMAT0027504	YP02119388
A12	hsa-let-7c-3p	MIMAT0026472	YP00205202	I12	hsa-miR-6803-3p	MIMAT0027507	YP02115732
A13	hsa-miR-1180-5p	MIMAT0026735	YP02100476	I13	hsa-miR-6804-3p	MIMAT0027509	YP02108252
A14	hsa-miR-1199-3p	MIMAT0031120	YP02102766	I14	hsa-miR-6804-5p	MIMAT0027508	YP02119084
A15	hsa-miR-1250-3p	MIMAT0026740	YP02110614	I15	hsa-miR-6805-5p	MIMAT0027510	YP02112777
A16	hsa-miR-1251-3p	MIMAT0026741	YP02119283	I16	hsa-miR-6806-3p	MIMAT0027513	YP02110431
A17	hsa-miR-1252-3p	MIMAT0026744	YP02114253	I17	hsa-miR-6806-5p	MIMAT0027512	YP02115030
A18	hsa-miR-1266-3p	MIMAT0026742	YP02109017	I18	hsa-miR-6807-3p	MIMAT0027515	YP02105778
A19	hsa-miR-1273h-3p	MIMAT0030416	YP02111365	I19	hsa-miR-6807-5p	MIMAT0027514	YP02119459
A20	hsa-miR-128-1-5p	MIMAT0026477	YP02117210	I20	hsa-miR-6808-3p	MIMAT0027517	YP02118534
A21	hsa-miR-128-2-5p	MIMAT0031095	YP02110351	I21	hsa-miR-6808-5p	MIMAT0027516	YP02106511
A22	hsa-miR-1288-5p	MIMAT0026743	YP02110095	I22	hsa-miR-6809-3p	MIMAT0027519	YP02111411
A23	hsa-miR-1296-3p	MIMAT0026637	YP02102262	I23	hsa-miR-6809-5p	MIMAT0027518	YP02101251
A24	hsa-miR-1298-3p	MIMAT0026641	YP02108006	I24	hsa-miR-6810-5p	MIMAT0027520	YP02101678
B01	hsa-miR-1301-5p	MIMAT0026639	YP02118512	J01	hsa-miR-6811-3p	MIMAT0027523	YP02111254
B02	hsa-miR-133a-5p	MIMAT0026478	YP00205501	J02	hsa-miR-6811-5p	MIMAT0027522	YP02114988
B03	hsa-miR-1343-5p	MIMAT0027038	YP02100503	J03	hsa-miR-6812-5p	MIMAT0027524	YP02110112
B04	hsa-miR-134-3p	MIMAT0026481	YP02114563	J04	hsa-miR-6813-5p	MIMAT0027526	YP02113477
B05	hsa-miR-1468-3p	MIMAT0026638	YP02110580	J05	hsa-miR-6814-5p	MIMAT0027528	YP02105238
B06	hsa-miR-152-5p	MIMAT0026479	YP02108962	J06	hsa-miR-6815-3p	MIMAT0027531	YP02102659
B07	hsa-miR-1537-5p	MIMAT0026765	YP02116182	J07	hsa-miR-6815-5p	MIMAT0027530	YP02106133
B08	hsa-miR-181d-3p	MIMAT0026608	YP02105350	J08	hsa-miR-6816-3p	MIMAT0027533	YP02109792
B09	gga-miR-190a-3p	MIMAT0026525	YP02100791	J09	hsa-miR-6816-5p	MIMAT0027532	YP02118525
B10	hsa-miR-1910-3p	MIMAT0026917	YP02110313	J10	hsa-miR-6817-5p	MIMAT0027534	YP02119229
B11	hsa-miR-208a-5p	MIMAT0026474	YP00205398	J11	hsa-miR-6818-3p	MIMAT0027537	YP02116125
B12	hsa-miR-208b-5p	MIMAT0026722	YP02110745	J12	hsa-miR-6818-5p	MIMAT0027536	YP02111698
B13	hsa-miR-210-5p	MIMAT0026475	YP02104321	J13	hsa-miR-6819-5p	MIMAT0027538	YP02110218
B14	hsa-miR-215-3p	MIMAT0026476	YP02114252	J14	hsa-miR-6820-3p	MIMAT0027541	YP02101831
B15	hsa-miR-216b-3p	MIMAT0026721	YP02117492	J15	hsa-miR-6820-5p	MIMAT0027540	YP02116855
B16	hsa-miR-2276-5p	MIMAT0026921	YP02108484	J16	hsa-miR-6821-3p	MIMAT0027543	YP02115191
B17	hsa-miR-3151-3p	MIMAT0027026	YP02118703	J17	hsa-miR-6821-5p	MIMAT0027542	YP02108669
B18	hsa-miR-3192-3p	MIMAT0027027	YP02107026	J18	hsa-miR-6822-3p	MIMAT0027545	YP02107987
B19	hsa-miR-328-5p	MIMAT0026486	YP02113205	J19	hsa-miR-6822-5p	MIMAT0027544	YP02112115
B20	hsa-miR-329-5p	MIMAT0026555	YP02107092	J20	hsa-miR-6823-3p	MIMAT0027547	YP02118078
B21	hsa-miR-370-5p	MIMAT0026483	YP02107400	J21	hsa-miR-6823-5p	MIMAT0027546	YP02114286
B22	hsa-miR-372-5p	MIMAT0026484	YP02115668	J22	hsa-miR-6824-3p	MIMAT0027549	YP02111157
B23	hsa-miR-383-3p	MIMAT0026485	YP02108841	J23	hsa-miR-6824-5p	MIMAT0027548	YP02102777
B24	hsa-miR-3912-5p	MIMAT0027036	YP02119128	J24	hsa-miR-6825-3p	MIMAT0027551	YP02115710
C01	hsa-miR-3928-5p	MIMAT0027037	YP02116032	K01	hsa-miR-6825-5p	MIMAT0027550	YP02104934

Well	miRNA ID	Accession #	Assay Catalog #	Well	miRNA ID	Accession #	Assay Catalog #
C02	hsa-miR-410-5p	MIMAT0026558	YP02111083	K02	hsa-miR-6826-3p	MIMAT0027553	YP02109244
C03	hsa-miR-412-5p	MIMAT0026557	YP02115902	K03	hsa-miR-6826-5p	MIMAT0027552	YP02102567
C04	hsa-miR-433-5p	MIMAT0026554	YP00205430	K04	hsa-miR-6827-3p	MIMAT0027555	YP02118217
C05	hsa-miR-4433b-3p	MIMAT0030414	YP02115942	K05	hsa-miR-6827-5p	MIMAT0027554	YP02114959
C06	hsa-miR-4433b-5p	MIMAT0030413	YP02118899	K06	hsa-miR-6828-3p	MIMAT0027557	YP02117525
C07	hsa-miR-487a-5p	MIMAT0026559	YP02112873	K07	hsa-miR-6828-5p	MIMAT0027556	YP02101120
C08	hsa-miR-487b-5p	MIMAT0026614	YP02115384	K08	hsa-miR-6829-3p	MIMAT0027559	YP02109974
C09	hsa-miR-489-5p	MIMAT0026605	YP02107863	K09	hsa-miR-6829-5p	MIMAT0027558	YP02107920
C10	hsa-miR-494-5p	MIMAT0026607	YP02111151	K10	hsa-miR-6830-3p	MIMAT0027561	YP02108622
C11	hsa-miR-500b-3p	MIMAT0027032	YP02105743	K11	hsa-miR-6830-5p	MIMAT0027560	YP02116880
C12	hsa-miR-6889-3p	MIMAT0027679	YP02104407	K12	hsa-miR-6831-3p	MIMAT0027563	YP02109795
C13	hsa-miR-511-3p	MIMAT0026606	YP02108582	K13	hsa-miR-6831-5p	MIMAT0027562	YP02109034
C14	hsa-miR-513b-3p	MIMAT0026749	YP02102666	K14	hsa-miR-6832-3p	MIMAT0027565	YP02116095
C15	hsa-miR-5189-3p	MIMAT0027088	YP02106506	K15	hsa-miR-6832-5p	MIMAT0027564	YP02112352
C16	hsa-miR-520g-5p	MIMAT0026611	YP02118483	K16	hsa-miR-6833-3p	MIMAT0027567	YP02119433
C17	hsa-miR-548ba	MIMAT0031175	YP02113326	K17	hsa-miR-6833-5p	MIMAT0027566	YP02114308
C18	hsa-miR-548e-5p	MIMAT0026736	YP02117828	K18	hsa-miR-6834-3p	MIMAT0027569	YP02107263
C19	hsa-miR-548f-5p	MIMAT0026739	YP02112172	K19	hsa-miR-6834-5p	MIMAT0027568	YP02100744
C20	hsa-miR-548j-3p	MIMAT0026737	YP02119385	K20	hsa-miR-6835-3p	MIMAT0027571	YP02116069
C21	hsa-miR-5699-5p	MIMAT0027103	YP02113295	K21	hsa-miR-6835-5p	MIMAT0027570	YP02119658
C22	hsa-miR-579-5p	MIMAT0026616	YP02117190	K22	hsa-miR-6836-3p	MIMAT0027575	YP02119660
C23	hsa-miR-580-5p	MIMAT0026617	YP02104524	K23	hsa-miR-6836-5p	MIMAT0027574	YP02105775
C24	hsa-miR-585-5p	MIMAT0026618	YP02114778	K24	hsa-miR-6837-3p	MIMAT0027577	YP02111870
D01	hsa-miR-598-5p	MIMAT0026620	YP02113869	L01	hsa-miR-6837-5p	MIMAT0027576	YP02118732
D02	hsa-miR-605-3p	MIMAT0026621	YP02105161	L02	hsa-miR-6838-3p	MIMAT0027579	YP02108844
D03	hsa-miR-619-5p	MIMAT0026622	YP02118098	L03	hsa-miR-6838-5p	MIMAT0027578	YP02109467
D04	hsa-miR-651-3p	MIMAT0026624	YP02112173	L04	hsa-miR-6839-3p	MIMAT0027581	YP02100700
D05	hsa-miR-6516-3p	MIMAT0030418	YP02101857	L05	hsa-miR-6839-5p	MIMAT0027580	YP02118949
D06	hsa-miR-6516-5p	MIMAT0030417	YP02117973	L06	hsa-miR-6840-3p	MIMAT0027583	YP02101589
D07	hsa-miR-653-3p	MIMAT0026625	YP02102836	L07	hsa-miR-6840-5p	MIMAT0027582	YP02110643
D08	hsa-miR-655-5p	MIMAT0026626	YP02104256	L08	hsa-miR-6841-3p	MIMAT0027585	YP02118210
D09	hsa-miR-656-5p	MIMAT0026627	YP02101477	L09	hsa-miR-6841-5p	MIMAT0027584	YP02107704
D10	hsa-miR-668-5p	MIMAT0026636	YP02108164	L10	hsa-miR-6842-3p	MIMAT0027587	YP02115310
D11	hsa-miR-670-3p	MIMAT0026640	YP02118565	L11	hsa-miR-6842-5p	MIMAT0027586	YP02118059
D12	hsa-miR-6889-5p	MIMAT0027678	YP02102772	L12	hsa-miR-6843-3p	MIMAT0027588	YP02105311
D13	hsa-miR-6726-3p	MIMAT0027354	YP02105180	L13	hsa-miR-6890-3p	MIMAT0027681	YP02102771
D14	hsa-miR-6726-5p	MIMAT0027353	YP02117526	L14	hsa-miR-6845-5p	MIMAT0027590	YP02110607
D15	hsa-miR-6728-3p	MIMAT0027358	YP02114170	L15	hsa-miR-6846-3p	MIMAT0027593	YP02107207
D16	hsa-miR-6730-5p	MIMAT0027361	YP02108676	L16	hsa-miR-6846-5p	MIMAT0027592	YP02106013
D17	hsa-miR-6731-3p	MIMAT0027364	YP02119123	L17	hsa-miR-6847-3p	MIMAT0027595	YP02103437
D18	hsa-miR-6732-3p	MIMAT0027366	YP02114748	L18	hsa-miR-6847-5p	MIMAT0027594	YP02113204
D19	hsa-miR-6732-5p	MIMAT0027365	YP02119428	L19	hsa-miR-6848-3p	MIMAT0027597	YP02101654
D20	hsa-miR-6733-3p	MIMAT0027368	YP02115361	L20	hsa-miR-6848-5p	MIMAT0027596	YP02107929
D21	hsa-miR-6733-5p	MIMAT0027367	YP02110920	L21	hsa-miR-6849-3p	MIMAT0027599	YP02109972
D22	hsa-miR-6734-3p	MIMAT0027370	YP02108135	L22	hsa-miR-6849-5p	MIMAT0027598	YP02116737
D23	hsa-miR-6734-5p	MIMAT0027369	YP02117038	L23	hsa-miR-6850-3p	MIMAT0027601	YP02113302
D24	hsa-miR-6735-5p	MIMAT0027371	YP02109568	L24	hsa-miR-6851-3p	MIMAT0027603	YP02117632
E01	hsa-miR-6736-3p	MIMAT0027374	YP02117652	M01	hsa-miR-6851-5p	MIMAT0027602	YP02109092
E02	hsa-miR-6736-5p	MIMAT0027373	YP02107530	M02	hsa-miR-6852-3p	MIMAT0027605	YP02116751
E03	hsa-miR-6737-5p	MIMAT0027375	YP02103653	M03	hsa-miR-6852-5p	MIMAT0027604	YP02104799

Well	miRNA ID	Accession #	Assay Catalog #	Well	miRNA ID	Accession #	Assay Catalog #
E04	hsa-miR-6738-3p	MIMAT0027378	YP02101750	M04	hsa-miR-6853-3p	MIMAT0027607	YP02118391
E05	hsa-miR-6738-5p	MIMAT0027377	YP02117886	M05	hsa-miR-6853-5p	MIMAT0027606	YP02105449
E06	hsa-miR-6739-3p	MIMAT0027380	YP02102392	M06	hsa-miR-6854-3p	MIMAT0027609	YP02108291
E07	hsa-miR-6739-5p	MIMAT0027379	YP02104637	M07	hsa-miR-6854-5p	MIMAT0027608	YP02115043
E08	hsa-miR-6740-3p	MIMAT0027382	YP02115616	M08	hsa-miR-6855-3p	MIMAT0027611	YP02112157
E09	hsa-miR-6740-5p	MIMAT0027381	YP02111295	M09	hsa-miR-6855-5p	MIMAT0027610	YP02116164
E10	hsa-miR-6741-3p	MIMAT0027384	YP02119233	M10	hsa-miR-6856-3p	MIMAT0027613	YP02113125
E11	hsa-miR-6741-5p	MIMAT0027383	YP02108759	M11	hsa-miR-6857-3p	MIMAT0027615	YP02116660
E12	hsa-miR-6742-3p	MIMAT0027386	YP02110516	M12	hsa-miR-6857-5p	MIMAT0027614	YP02111562
E13	hsa-miR-6742-5p	MIMAT0027385	YP02119213	M13	hsa-miR-6858-5p	MIMAT0027616	YP02112076
E14	hsa-miR-6743-5p	MIMAT0027387	YP02117389	M14	hsa-miR-6859-5p	MIMAT0027618	YP02103332
E15	hsa-miR-6744-3p	MIMAT0027390	YP02114950	M15	hsa-miR-6860	MIMAT0027622	YP02105127
E16	hsa-miR-6744-5p	MIMAT0027389	YP02104363	M16	hsa-miR-6861-5p	MIMAT0027623	YP02119257
E17	hsa-miR-6746-3p	MIMAT0027393	YP02116628	M17	hsa-miR-6862-3p	MIMAT0027626	YP02116509
E18	hsa-miR-6747-3p	MIMAT0027395	YP02113949	M18	hsa-miR-6862-5p	MIMAT0027625	YP02116492
E19	hsa-miR-6748-3p	MIMAT0027397	YP02106726	M19	hsa-miR-6863	MIMAT0027627	YP02112191
E20	hsa-miR-6748-5p	MIMAT0027396	YP02103103	M20	hsa-miR-6864-3p	MIMAT0027629	YP02105403
E21	hsa-miR-6749-5p	MIMAT0027398	YP02108544	M21	hsa-miR-6864-5p	MIMAT0027628	YP02106668
E22	hsa-miR-6750-5p	MIMAT0027400	YP02111905	M22	hsa-miR-6865-3p	MIMAT0027631	YP02115783
E23	hsa-miR-6751-3p	MIMAT0027403	YP02106215	M23	hsa-miR-6865-5p	MIMAT0027630	YP02118138
E24	hsa-miR-6751-5p	MIMAT0027402	YP02118769	M24	hsa-miR-6866-3p	MIMAT0027633	YP02118145
F01	hsa-miR-6752-5p	MIMAT0027404	YP02110467	N01	hsa-miR-6866-5p	MIMAT0027632	YP02111715
F02	hsa-miR-6753-3p	MIMAT0027407	YP02118515	N02	hsa-miR-6867-3p	MIMAT0027635	YP02111051
F03	hsa-miR-6753-5p	MIMAT0027406	YP02105614	N03	hsa-miR-6867-5p	MIMAT0027634	YP02116720
F04	hsa-miR-6754-5p	MIMAT0027408	YP02116950	N04	hsa-miR-6868-3p	MIMAT0027637	YP02102933
F05	hsa-miR-6755-3p	MIMAT0027411	YP02117866	N05	hsa-miR-6868-5p	MIMAT0027636	YP02111215
F06	hsa-miR-6755-5p	MIMAT0027410	YP02109049	N06	hsa-miR-6869-3p	MIMAT0027639	YP02103497
F07	hsa-miR-6756-5p	MIMAT0027412	YP02119429	N07	hsa-miR-6870-3p	MIMAT0027641	YP02117175
F08	hsa-miR-6757-3p	MIMAT0027415	YP02107562	N08	hsa-miR-6871-3p	MIMAT0027643	YP02103093
F09	hsa-miR-6757-5p	MIMAT0027414	YP02109195	N09	hsa-miR-6871-5p	MIMAT0027642	YP02113232
F10	hsa-miR-6758-3p	MIMAT0027417	YP02109503	N10	hsa-miR-6872-5p	MIMAT0027644	YP02101577
F11	hsa-miR-6759-3p	MIMAT0027419	YP02115722	N11	hsa-miR-6873-3p	MIMAT0027647	YP02102681
F12	hsa-miR-6759-5p	MIMAT0027418	YP02107923	N12	hsa-miR-6873-5p	MIMAT0027646	YP02103510
F13	hsa-miR-6761-3p	MIMAT0027423	YP02106554	N13	hsa-miR-6874-3p	MIMAT0027649	YP02117460
F14	hsa-miR-6762-3p	MIMAT0027425	YP02106317	N14	hsa-miR-6874-5p	MIMAT0027648	YP02109241
F15	hsa-miR-6762-5p	MIMAT0027424	YP02105421	N15	hsa-miR-6875-3p	MIMAT0027651	YP02108527
F16	hsa-miR-6763-5p	MIMAT0027426	YP02118635	N16	hsa-miR-6875-5p	MIMAT0027650	YP02116574
F17	hsa-miR-6764-3p	MIMAT0027429	YP02117884	N17	hsa-miR-6876-3p	MIMAT0027653	YP02115467
F18	hsa-miR-6764-5p	MIMAT0027428	YP02109894	N18	hsa-miR-6876-5p	MIMAT0027652	YP02110823
F19	hsa-miR-6765-5p	MIMAT0027430	YP02114389	N19	hsa-miR-6877-5p	MIMAT0027654	YP02115155
F20	hsa-miR-6766-5p	MIMAT0027432	YP02102746	N20	hsa-miR-6878-3p	MIMAT0027657	YP02110113
F21	hsa-miR-6767-3p	MIMAT0027435	YP02109520	N21	hsa-miR-6878-5p	MIMAT0027656	YP02113438
F22	hsa-miR-6767-5p	MIMAT0027434	YP02118572	N22	hsa-miR-6879-5p	MIMAT0027658	YP02114612
F23	hsa-miR-6768-3p	MIMAT0027437	YP02102014	N23	hsa-miR-6880-5p	MIMAT0027660	YP02116477
F24	hsa-miR-6768-5p	MIMAT0027436	YP02110713	N24	hsa-miR-6881-3p	MIMAT0027663	YP02101083
G01	hsa-miR-6769a-5p	MIMAT0027438	YP02106693	O01	hsa-miR-6881-5p	MIMAT0027662	YP02109019
G02	hsa-miR-6769b-3p	MIMAT0027621	YP02101670	O02	hsa-miR-6882-3p	MIMAT0027665	YP02108189
G03	hsa-miR-6769b-5p	MIMAT0027620	YP02113037	O03	hsa-miR-6882-5p	MIMAT0027664	YP02106235
G04	hsa-miR-6770-3p	MIMAT0027441	YP02111845	O04	hsa-miR-6883-3p	MIMAT0027667	YP02117395
G05	hsa-miR-6770-5p	MIMAT0027440	YP02103217	O05	hsa-miR-6883-5p	MIMAT0027666	YP02118150

Well	miRNA ID	Accession #	Assay Catalog #	Well	miRNA ID	Accession #	Assay Catalog #
G06	hsa-miR-6771-3p	MIMAT0027443	YP02119355	O06	hsa-miR-6884-3p	MIMAT0027669	YP02107360
G07	hsa-miR-6771-5p	MIMAT0027442	YP02105212	O07	hsa-miR-6884-5p	MIMAT0027668	YP02118648
G08	hsa-miR-6772-3p	MIMAT0027445	YP02103982	O08	hsa-miR-6885-3p	MIMAT0027671	YP02101345
G09	hsa-miR-6772-5p	MIMAT0027444	YP02118541	O09	hsa-miR-6885-5p	MIMAT0027670	YP02119482
G10	hsa-miR-6773-3p	MIMAT0027447	YP02106901	O10	hsa-miR-6886-5p	MIMAT0027672	YP02105593
G11	hsa-miR-6773-5p	MIMAT0027446	YP02116984	O11	hsa-miR-6887-3p	MIMAT0027675	YP02116052
G12	hsa-miR-6774-3p	MIMAT0027449	YP02106422	O12	hsa-miR-6887-5p	MIMAT0027674	YP02113868
G13	hsa-miR-6774-5p	MIMAT0027448	YP02107421	O13	hsa-miR-6888-3p	MIMAT0027677	YP02111557
G14	hsa-miR-6775-3p	MIMAT0027451	YP02118761	O14	hsa-miR-6888-5p	MIMAT0027676	YP02104718
G15	hsa-miR-6775-5p	MIMAT0027450	YP02119417	O15	hsa-miR-6890-5p	MIMAT0027680	YP02110989
G16	hsa-miR-6776-3p	MIMAT0027453	YP02114654	O16	hsa-miR-6891-3p	MIMAT0027683	YP02106363
G17	hsa-miR-6776-5p	MIMAT0027452	YP02116812	O17	hsa-miR-6891-5p	MIMAT0027682	YP02102735
G18	hsa-miR-6777-5p	MIMAT0027454	YP02101311	O18	hsa-miR-6892-3p	MIMAT0027685	YP02114726
G19	hsa-miR-6778-3p	MIMAT0027457	YP02105003	O19	hsa-miR-6892-5p	MIMAT0027684	YP02113175
G20	hsa-miR-6778-5p	MIMAT0027456	YP02101666	O20	hsa-miR-6893-3p	MIMAT0027687	YP02119432
G21	hsa-miR-6779-3p	MIMAT0027459	YP02114622	O21	hsa-miR-6893-5p	MIMAT0027686	YP02116619
G22	hsa-miR-6779-5p	MIMAT0027458	YP02117843	O22	hsa-miR-7107-3p	MIMAT0028112	YP02101788
G23	hsa-miR-6780a-3p	MIMAT0027461	YP02117483	O23	hsa-miR-7108-3p	MIMAT0028114	YP02101501
G24	hsa-miR-6780a-5p	MIMAT0027460	YP02117364	O24	hsa-miR-7109-5p	MIMAT0028115	YP02111712
H01	hsa-miR-6780b-3p	MIMAT0027573	YP02100688	P01	hsa-miR-7110-3p	MIMAT0028118	YP02114244
H02	hsa-miR-6781-5p	MIMAT0027462	YP02116752	P02	hsa-miR-7111-3p	MIMAT0028120	YP02110682
H03	hsa-miR-6782-3p	MIMAT0027465	YP02104981	P03	hsa-miR-7112-3p	MIMAT0028122	YP02106857
H04	hsa-miR-6783-3p	MIMAT0027467	YP02111036	P04	hsa-miR-7112-5p	MIMAT0028121	YP02109316
H05	hsa-miR-6783-5p	MIMAT0027466	YP02113087	P05	hsa-miR-7113-3p	MIMAT0028124	YP02106063
H06	hsa-miR-6785-3p	MIMAT0027471	YP02118992	P06	hsa-miR-7113-5p	MIMAT0028123	YP02107502
H07	hsa-miR-6785-5p	MIMAT0027470	YP02103050	P07	hsa-miR-7114-3p	MIMAT0028126	YP02109524
H08	hsa-miR-6786-3p	MIMAT0027473	YP02115225	P08	hsa-miR-7114-5p	MIMAT0028125	YP02115348
H09	hsa-miR-6786-5p	MIMAT0027472	YP02107023	P09	hsa-miR-7150	MIMAT0028211	YP02109509
H10	hsa-miR-6787-3p	MIMAT0027475	YP02104871	P10	hsa-miR-7151-3p	MIMAT0028213	YP02108098
H11	hsa-miR-6787-5p	MIMAT0027474	YP02119387	P11	hsa-miR-7151-5p	MIMAT0028212	YP02113505
H12	hsa-miR-6788-3p	MIMAT0027477	YP02106685	P12	BlankSpot	N/A	YP02119291
H13	hsa-miR-6788-5p	MIMAT0027476	YP02107621	P13	SNORD44 (hsa)	N/A	YP00203902
H14	hsa-miR-6789-5p	MIMAT0027478	YP02100084	P14	SNORD38B (hsa)	N/A	YP00203901
H15	hsa-miR-6790-3p	MIMAT0027481	YP02105201	P15	SNORD49A (hsa)	N/A	YP00203904
H16	hsa-miR-6790-5p	MIMAT0027480	YP02105097	P16	U6 snRNA (v2)	N/A	YP02119464
H17	hsa-miR-6791-3p	MIMAT0027483	YP02113917	P17	cel-miR-39-3p	MIMAT0000010	YP00203952
H18	hsa-miR-6792-3p	MIMAT0027485	YP02115808	V	UniSp2	N/A	YP00203950
H19	hsa-miR-6792-5p	MIMAT0027484	YP02117927	P19	UniSp4	N/A	YP00203953
H20	hsa-miR-6793-5p	MIMAT0027486	YP02114368	P20	UniSp5	N/A	YP00203955
H21	hsa-miR-6794-3p	MIMAT0027489	YP02102949	P21	UniSp6	N/A	YP00203954
H22	hsa-miR-6794-5p	MIMAT0027488	YP02116385	P22	UniSP3	N/A	YP02119288
H23	hsa-miR-6795-3p	MIMAT0027491	YP02119119	P23	UniSP3	N/A	YP02119288
H24	hsa-miR-6795-5p	MIMAT0027490	YP02103188	P24	UniSP3	N/A	YP02119288

Ordering Information

Product	Contents	Cat. no.
miRCURY LNA miRNA Focus PCR Panels	miRCURY LNA miRNA PCR Panels for application-based miRNome profiling, available in 96-well or 384-well format; for SYBR® Green-based detection	339325
miRCURY LNA miRNA miRNome PCR Panels	miRCURY LNA miRNA PCR Panels for PCR-based miRNome profiling, available in 384-well format; for SYBR® Green-based detection	339322
miRCURY LNA miRNA QC PCR Panel	miRCURY LNA miRNA PCR Panel of quality control assays, available in 96-well or 384-well format; for SYBR® Green-based detection	339331
miRCURY LNA miRNA Custom PCR Panels	8 identical, ready-to-use 96- or 384-well plates; each well contains primers sufficient for one 10 µl reaction; for SYBR® Green-based detection	339330
miRCURY LNA Custom PCR Panel Additional Plate	Additional miRCURY LNA Custom PCR Panel plates; set of 4 plates; only available in addition to the base plates ordered through the core product (cat. no. 339330)	339332
miRCURY LNA miRNA PCR Assays	Contains forward and reverse primers for 200 SYBR® Green-based, real-time qPCR reactions, 166 EvaGreen-based digital PCR reactions for Nanoplate 8.5k or 50 EvaGreen-based digital PCR reactions for Nanoplate 26k	339306
miRCURY LNA miRNA Probe PCR Assays	Complete premixed assays containing LNA-enhanced target-specific forward primer and probe. For 200 reactions.	339350
miRCURY LNA miRNA Custom Probe PCR Assays	Custom-designed, target-specific forward primer and probe for any user-defined miRNA target. Complete premixed assay for 200 reactions.	339351

Related Products

Product	Contents	Cat. no.
miRCURY LNA RT Kit	For 8–64 cDNA synthesis reactions: 5x RT SYBR Green Reaction Buffer, 5x RT Probe Reaction Buffer, 10x RT Enzyme Mix, UniSp6, RNA Spike-in template, RNase-Free Water	339340
RNA Spike-In Kit, For RT	Contains the UniSp2, UniSp4, and UniSp5 RNA Spike-in Template Mix and the cel-miR-39-3p RNA Spike-in Template	339390
miRCURY LNA SYBR® Green PCR Kits (200)	For 200 reactions: 2X miRCURY SYBR Green Master Mix, RNase-Free Water	339345
miRCURY LNA SYBR® Green PCR Kits (600)	For 600 reactions: 2X miRCURY SYBR Green Master Mix, RNase-Free Water	339346
miRCURY LNA SYBR® Green PCR Kits (4000)	For 4000 reactions: 2X miRCURY SYBR Green Master Mix, RNase-Free Water	339347
miRCURY LNA Probe PCR Kit (200)	For 200 reactions: 2X QuantiNova Probe Master Mix, 10X miRCURY Probe Univ. Primer, Rox Reference Dye, RNase-Free Water	339371
miRCURY LNA Probe PCR Kit (800)	For 800 reactions: 2X QuantiNova Probe Master Mix, 10X miRCURY Probe Univ. Primer, Rox Reference Dye, RNase-Free Water	339372
miRCURY LNA Probe PCR Kit (4000)	For 4000 reactions: 2X QuantiNova Probe Master Mix, 10X miRCURY Probe Univ. Primer, Rox Reference Dye, RNase-Free Water	339373
miRCURY LNA miRNA PCR Starter Kit	Two miRCURY LNA PCR Assays of your choice, spike-in control Assay (UniSp6), one candidate endogenous control assay (miR-103-3p) and all reagents for 20 reverse transcription reactions and 100 PCR amplifications; for SYBR® Green-based qPCR detection	339320

Important

The miRCURY LNA™ miRNA PCR Panels are Ready-to-Use and designed for optimal performance with the miRCURY LNA RT Kit and the miRCURY LNA SYBR® Green PCR Kit. The performance of the primer sets will be affected when used in combination with less than optimal reagents. Use the miRCURY LNA miRNA PCR Panel Handbook for experiment setup. RNA work requires specific handling and precautions should be taken to prevent RNase contamination and degradation of the RNA sample and reagents.

Shipping and storage

The Ready-to-Use miRCURY LNA miRNA PCR Panels are shipped at room temperature and can be stored at 4°C for at least 6 months. For long term storage, it is recommended to place the panels at -20°C. Under these conditions, the LNA PCR primers are stable for at least 6 months after receipt.

Intended use

miRCURY LNA miRNA PCR Assays are intended for molecular biology applications. This product is not intended for the diagnosis, prevention or treatment of a disease.

Safety information

When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, please consult the appropriate safety data sheets (SDSs). These are available online in convenient and compact PDF format at www.qiagen.com/safety where you can find, view and print the SDS for each QIAGEN kit and kit component.

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual. QIAGEN kit handbooks and user manuals are available at www.qiagen.com or can be requested from QIAGEN Technical Services or your local distributor.

Trademarks: QIAGEN®, Sample to Insight®, GeneGlobe®, LNA™ (QIAGEN Group), SYBR® (Life Technologies Corporation). Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law. 1107212 05/2017 HB-2376-001 © 2017 QIAGEN, all rights reserved.
